



Ramaciotti Centre for Genomics

Guide to Online Submission of Samples

LIMS Submission Guide

Ramaciotti Centre for Genomics

Biosciences South (E26) UNSW Sydney Kensington NSW 2052 Australia

ramaciotti.unsw.edu.au (02) 9385 1241 ramaciotti@unsw.edu.au



Australia's Global University

- 1. Navigate to the sample submission portal: <u>https://ramaciotti.clarityLIMS.com/lablink/Welcome.do</u>
- 2. Log onto LIMS using your User ID and password. New users must register for a User ID.
- 3. Click on Get Sample Submission Forms and Instructions.



4. Choose the appropriate form for the service you require and download.

R		Your Profile	Change Your Pasaword	Request a User ID		Sign.Qut
Remucistti Centre for Genomics	Download Files		System Admi Jak 31, 3018	nistratur at Administrative L	ab ide not pickl sig	ped in as admin
LubLink Home					Get help f	or this page
Submit Samples	Example Sample Sheet, Example only. Please do not use Biomarkger, Sample Submission Form Illiomad, Sample Submission Form					
Get Samele Submasia Farms and Instructions	Initiania kuka: Untegela Auditata Initiania kuka: Untegela Auditata Initiania kuka: Untegela Auditata Initiania kuka: Untegela Auditatata Initiania kuka: Untegela Auditatatatatatatatatatatatatatatatatatata					
Open the Lablick Hela	Eludigen Samtle Submission Form WanoString Samtle Submission Form					

- **5.** Fill in <u>all sections</u> of the form paying close attention to the instructions located at the beginning of the form.
- 6. Save the form to an appropriate directory on your computer.

Important:

- You cannot submit samples unless you have a Purchase Order (PO) or a UNSW account code (internal UNSW users only). Your finance department can generate a PO against the quote we have issued for your project.
- Do not use duplicate sample names.
- Use only alphanumerical characters and underscores in the sample name.

7. Click on Submit Samples.

	R				Your Profile	Change Your Password	Request a User ID	Contact.Us	Sign.Out
		me Page - Project List					Youla Russell at Ran Jul 13, 2013	nacietti Centre sigr	ed in as Taniall
Click here to submit	LabLink Hume	Trejects Samples Files							
samples.	Submit Samples							C Get help for	this page
	Con Set Sample Submission	View prejects with open		Appry					
	1	Name	+ LIMSED		Status	Date Greated			
	Open the LabLink Help								
	100 m								
	Property by Canadatana								

8. Create a project name. Click Next.

Important:

- Please include your surname in the project name.
- Choose a project name that will help to differentiate the project from any past or future projects you have/plan to submit.

(K)				<u>Your Profile</u>	Change Your Password	<u>Request a User ID</u>	<u>Contact Us</u>	Sign Out
Ramaciotti Centre for Genomics	Sample Submission - Creat	e a New Project				Tonia Russell at Ran Jul 13, 2015	naciotti Centre sign	ed in as ToniaR
LabLink Home			Next	Cancel				
Submit Samples Get Sample Submission Forms and Instructions	Required fields have	a red outline.					@ Get help fo	or this page
Porms and instructions		Enter Project Information					-	
		Project name:						
Open the LabLink Help		Note:	·]	
							4	
			Next	Cance	A Committee			

9. Upload the completed submission form that you have saved on your computer. Click Next.

				Your Profile	Giange Tour Password	Request a User ID	Contact Us	Sign Out
Ramaciotti Centre for Genomics	Enter Sample Informatio	n for Project - ghyhjhy				Tonia Russell at Ran Jul 12, 2010	nacietti Centre sign	ed in as ToniaR
abLink Home			Next	Cancel				
ubmit Samples							() Get help fo	or this page
Get Sample Submission Forms and Instructions		Choose File No file chosen	C				Ē.	
		The maximum file size that						
Open the <u>LabLink Help</u>		L						
J tile			Next	Cancel				

10. You will be asked to review your sample details and agree to the Ramaciotti Centre Terms and Conditions (Figure A). After clicking agree, you will be prompted to print out your submission form (Figure B). Please print, sign and send to the Ramaciotti Centre with your samples. This form also contains the assigned LIMS sample IDs.

Review You	ur Samples												Tonia Ru Jul 13, 20	ssell at Ramaciotti Con	tre signed in as To
Please re	view your s	ample inf	ormation.	Click Agree t	o signify	your understa	nding and ad	ceptance o	of our Term	s and Condit	ions.				
										2010 121 20				@ Get !	help for this p
Namet	LIMS :	Status‡	Project¢	Date Received	Last Note	Application ¢	Progress¢	Sample : Buffer	Sample Conc. Ing per ull	<u>Platform</u> ≎	Tissue Type	Extraction : Method	Purchase Order_UNSW¢ ACC Code	Quote No. 🜩	Research Discipline
ABC	RUS549A1	Pending	ghyhjhy			20Kb Genomic		Tris HCL	250.00	PacBio	culture	Qiagen	12345	PacBioQ123JUL15	Biology (general)
DEF	RUS549A2	Pending	ghyhjhy			20Kb Genomic		Tris HCL	250.00	PacBio	culture	Giagen	12345	PacBio0123JUL15	Food/Nutritio Sciences
	Terms & Conditions of Service Use of the Ramaclotti Centre for Genomics (RCG) services and consumables are governed by the following conditions: Definitions: In these terms: (a) Data means the data or other results generated by RCG from the analysis of the Samples. (b) Cutote means the data or other results generated by RCG from the analysis of the Services and the pricing. (c) Samples means the samples provided by the institution to RCG Ord the data so the Services and the pricing. (d) Services means analysis and related services to be provided by RCG as indicated in the Submission Form and which is further described on the RCG website. (e) Submission Form means the document you completed and submitted setting out the services required, sample details and other relevant														
						_	Back	Disa	orma	Agree				·	

A – Agree to Terms & onditions

B – Print out form & sign

R										Y	our Profile	Change	Your Passwo	rd Request a l	Jser ID Con
ciotti Centre Genomica	Summary	of Submitte	ed Sample	IS .										Tonia Ru Jul 13, 20	sell at Ramaciotti (
ome	PLEASE	PRINT AND	SIGN THE	S PAGE. TH	IS FORM SH	OULD BE	E SENT WITH Y	OUR SAMPL	ES.						
ples	IMPORTA	NT:													
	Tube sub	missions m	ust be cle	arly labele	d with the be	low corn	esponding san	nple LIMS ID	(e.g. RUS	140A1). Do	not send you	r samples	unless they	are labeled with t	hese IDs.
ubmission structions	Plate sub these IDs		nould be l	abeled with	ı your contair	ner (plate	e) name (as sh	own on your	submissio	n form) an	d Project LIN	IS ID (e.g.	RUS140). Do	not send your pla	tes unless they
bLink Help															G G
ink Help	<u>Name</u> ¢	LIMS ID ¢	<u>Status</u> ¢	Project¢	Date Received	<u>Last</u> ¢ Note	<u>Application</u> ¢	Progress	<u>Sample</u> ¢ Buffer	Sample Conc. Ing per	<u>Platform</u> ≎	<u>Tissue</u> ¢ <u>Type</u>	Extraction Method	Purchase Order UNSW¢ ACC Code	<u>Quote No.</u>
	ABC	RUS549A1	Pending	ghyhjhy			20Kb Genomic		Tris HCL	250.00	PacBio	culture	Giagen	12345	PacBioQ123JUL
	DEE	RUS549A2	Pending	ghyhjhy			20Kb Genomic		Tris HCL	250.00	PacBio	culture	Giagen	12345	PacBioQ123JUL
	×	MS		ature		_									

- **11.** Please clearly label your tubes or plates with their corresponding LIMS sample ID (e.g. RUS549A1 in screen shot B above). Plates should also be labelled with their container ID e.g. RUS549 plate 1, RUS549 plate 2 etc.
- **12.** Once you have clicked the done button, the sample submission form has been submitted to the Centre.